## Bürkle GmbH

Rheinauen 5 | D - 79415 Bad Bellingen  
Telefon +49 (0) 76 35 8 27 95-0  
Telefax +49 (0) 76 35 8 27 95-31  
info@buerkle.de | www.buerkle.de

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## General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

## Safety instructions

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to escaping or off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood. Wear appropriate protective clothing.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ The instrument is suitable for indoor use and manual operation. Do not expose the device to any direct sunlight.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

### CAUTION!

Escaping liquids due to falling over or tipping of the instrument. Bodily injury due to escaping liquid and damage to the instrument.

- ▶ Make certain the container is positioned with adequate stability.
- ▶ Do not transport with the stopcock screwed on.

## Layout and function

- ▶ The device is made of PP, PE and PTFE.
- ▶ The device is suitable for dispensing non-crystallizing liquids (without solids) from canisters and barrels.
- ▶ The device can be used in the temperature range of 0 - 40 °C (briefly up to 60 °C).
- ▶ The device is pressure-resistant without a discharge up to max. 2 bar at 20 °C.

Point out any damage in the packaging to the freight forwarder/supplier immediately.

## Assembly and safe usage

### CAUTION!

Possible danger of breakage.

- ▶ Apply only normal manual force.
- ▶ Do not lift the canister by the stopcock.
- ▶ Do not suspend filling containers from the stopcock.

1. Insert the white PE seal ② in the sealing flange inside the union nut ①.
2. Screw the stop cock with union nut ① onto the container thread (R 3/4" connecting thread).
3. If applicable insert the discharge reducing nozzle ⑤.

### ! WARNING!

Danger of bodily injury due to liquid under high pressure escaping.

- ▶ The pressure in the canister must not exceed 2 bar.
- ▶ When the discharge reducing nozzle ⑤ is being used, only a non-pressurized discharge is permitted (nozzle is only inserted).

### CAUTION!

Openings and spaces on the device. Danger of crushing.

- ▶ Be careful for spaces and openings when activating the plug.

4. Open the plug ⑥ (turn anti-clockwise) and dispense liquid.
5. Close the plug (turn clockwise).
6. Turn the angle discharge ④ up after dispensing to prevent dripping. Do not tighten the union nut ③ too tightly when doing this.

### ! WARNING!

Bodily injury due to escaping liquid

- ▶ Close the stopcock completely after dispensing (set plug to at least „zero“).
- ▶ Keep the stopcock closed when it is not in use.
- ▶ Do not unscrew the stopcock from the container when it is filled.
- ▶ Empty the stopcock completely before disassembly.

## Troubleshooting

Problem	To eliminate the error
Thread connection is not sealed	<ol style="list-style-type: none"> <li>1. Insert PE seal.</li> <li>2. Thread is damaged (contact dealer or manufacturer).</li> </ol>
Stopcock is dripping	<ol style="list-style-type: none"> <li>1. Close the plug more tightly.</li> <li>2. Turn the discharge so it is pointing up.</li> </ol>

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