**Bürkle GmbH**

Rheinauen 5 | D - 79415 Bad Bellingen  
Telefon +49 (0) 76 35 8 27 95-0  
Telefax +49 (0) 76 35 8 27 95-31  
info@buerkle.de | www.buerkle.de

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### General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

### Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ Do not fill the instrument with potentially explosive substances.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.

#### **⚠ WARNING!**

Damage to the respiratory tract and other organs due to escaping media.

- ▶ When sampling hazardous substances, follow the safety data sheet of the medium to be sampled.
- ▶ Wear the appropriate protective clothing.

#### **NOTE!**

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
  - ▶ Pay attention to the material resistance.
- ▶ Do not use any force when operating the instrument.

### Layout and function

- ▶ The instrument is suitable for sampling small quantities of powders and granulates.
- ▶ The instrument is made of stainless steel 1.4404 (AISI 316 L) / 1.4571 (AISI 316Ti).
- ▶ The instrument has three sample chambers for one sample zone.

- ▶ Exchangeable sample inserts allow sample quantities from 0 up to 4.0 ml (see accessories).
- ▶ The instrument allows to withdraw a triple-sample without inserting the instrument repeatedly into the sampling material.
- ▶ The instrument is maintenance-free and suitable for steam sterilisation (up to 121°C).

### Parts in contact with medium

- ▶ Stainless steel 1.4404 (AISI 316 L) / 1.4571 (AISI 316Ti)

Point out any damage in the packaging to the freight forwarder/supplier immediately.

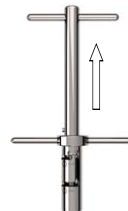
### Assembly

#### **CAUTION!**

Openings with sharp edges on the instrument. Danger of cutting.

- ▶ Keep hands well away from openings with sharp edges.

1. Set the instrument in the position for „disassembling“.

2. Pull the inner rod ④ of the instrument completely out.
3. Place the sample inserts into the inner rod ④. Sample 1 corresponds to „①“, Sample 2 corresponds to „②“ and Sample 3 corresponds to „③“.
4. Reinsert the inner rod ④ with the sample inserts into the sampling tube ⑤.

### Safe sampling

#### **CAUTION!**

Pointed parts on the instrument. Risk of injury.

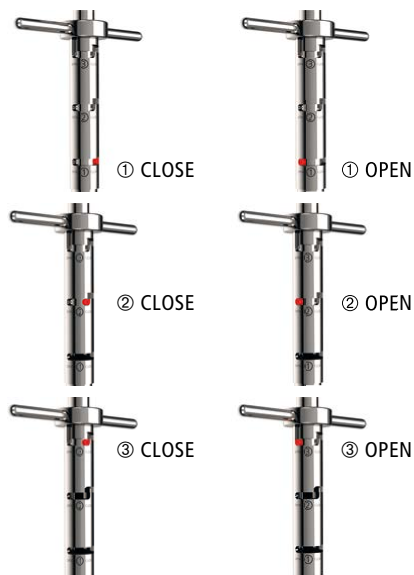
- ▶ Do not transport the instrument when the tip is not secured. Do not throw the instrument around.
- ▶ Do not pierce the instrument completely through the container.

#### **NOTE!**

The instrument is designed for manual operation.

- ▶ Do not hammer the instrument into the container by using a hammer.

1. Set the instrument at „①“ in position „CLOSE“.
2. The instrument is closed and can pierce into the powder/granulate.
3. Pierce the instrument into the powder/granulate up to the desired depth (the scale indicates the depth in cm).
4. Set the instrument at „①“ in position „OPEN“.
5. The instrument opens and the sample is filled.
6. Set the instrument at „①“ in position „CLOSE“.
7. The instrument is closed and can be removed from the powder/granulate.
8. If a second sample should be taken, set the instrument at „②“ in position „CLOSE“.
9. Set the instrument at „②“ in position „OPEN“. The instrument opens and is filled with the second sample.
10. Set the instrument at „②“ in position „CLOSE“.
11. The instrument is closed and can be removed from the powder/granulate.
12. If a third sample should be taken, set the instrument at „③“ in position „CLOSE“.
13. Set the instrument at „③“ in position „OPEN“. The instrument opens and is filled with the third sample.
14. Set the instrument at „③“ in position „CLOSE“.
15. The instrument is closed and can be removed from the powder/granulate.
16. Open the instrument and empty the sample on an appropriate surface or into a suitable collecting container.

#### After taking the sample

1. Set the instrument in the position for „disassembling“.
2. Pull the Inner rod ④ of the instrument completely out and remove the sample inserts.
3. Unscrew the tip ⑥.
4. Clean all parts. Use the cleaning brushes (see accessories) for the inside of tubes.

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